

# AMYLASE PAPER SALIVA

## AMYLASE MAPPING

- + fast and reliable presumptive test for saliva
- + on any type of fabric or material
- + compatible with other detection techniques

1. Moisten the Amylase Paper with distilled water
2. Place over the suspected stain and press for 10-15 seconds
3. Remove the Amylase Paper from the suspected stain and incubate for 10 minutes at room temperature
4. Pour or spray the Iodine working solution over the Amylase Paper

## INTERPRETATION OF RESULTS

**POSITIVE** a white area surrounded by a blue background indicates a presumptive positive result for saliva

**NEGATIVE** no white area is visible on the Amylase Paper

Presumptive results can be confirmed with an alpha-Amylase specific SERATEC® AmylaseTest: Simply collect/cut the area of the suspected stain, extract in the provided dilution buffer and run a SERATEC® AmylaseTest for more confidence.

**SERATEC®**  
EXCELLENCE IN FORENSIC BIOLOGY

SERATEC® GmbH  
P.O. Box 37 06  
37027 Goettingen  
Germany

Phone: +49 (0)551 504800  
Fax: +49 (0)551 5048080  
E-mail: [contact@seratec.com](mailto:contact@seratec.com)  
Web: [www.seratec.com](http://www.seratec.com)

